


Analysis of Tryptophan metabolites

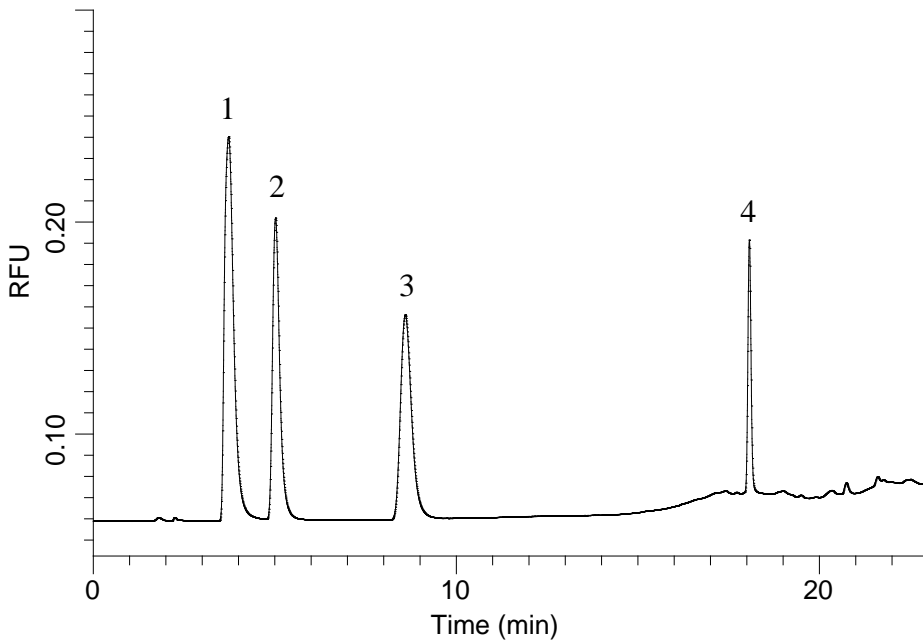

Conditions

System : Accustream system + LIF726
Column : Monocap C18 Fast-flow (150 x 0.2 mm I.D.)
Column Cat. No. : 5020-10111
Eluent : A) 10mM CH₃COONa (pH4.0, CH₃COOH)/CH₃OH = 99/1, v/v
 B) 10mM CH₃COONa (pH4.0, CH₃COOH)/CH₃OH = 7/93, v/v
 A/B = 99/1 - 6min - 83/17 - 2min - 30/70 - 12min - 30/70 - 0.1min - 99/1 - 5min - 99/1, v/v
Flow Rate : 1 μL/min
Col. Temp. : 30 °C
Detection : FL Ex 266 nm Em 300-350 nm (LIF726)
Injection Vol. : 0.02 μL
Sample : Standard

Analyte:

1. Serotonin	2.3mg/L
2. Tryptophan	1.9mg/L
3. 5-Hydroxyindoleacetic acid	2.4mg/L
4. Melatonin	1.6mg/L